4-Channel Timer

Cat. # 86-TMR-4

- Four independent countdown/count-up timers (T1–T4) with clock and alarm functions can run simultaneously; each has its own flashing indicator.
- Automatic count-up begins when a countdown reaches zero.
- Includes clock mode, 12/24-hour format, memory recall, and 1-minute buzzer alarm.

Clock Mode

- Enter Clock Mode: Press CLOCK. Time and flashing colon will display.
- Toggle 12/24hr Format: Hold START/STOP for 3 seconds.

Set Clock Time:

- Hold CLOCK for 3 seconds (until beep).
- "HOUR", "MINUTE", "SECOND" flash on display. "P" appears for PM in 12hr mode.
- Adjust using HOUR, MINUTE, SECOND (hold to fast set).
- Press the SECOND button to reset seconds:
 - If 00–29: resets to 00.
 - If 30–59: resets to 00 and adds 1 minute.
- Press CLOCK again to confirm and exit.

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Multi-Timer Function

- All four timers (T1–T4) can run simultaneously.
- While running, each timer's indicator flashes.
- At completion, the buzzer will sound and, the corresponding indicator flashes slower.

Countdown Timer (T1–T4)

- Press T1, T2, T3, or T4 to select a timer. Indicator appears on display
- Set time using HOUR, MINUTE, SECOND (hold for fast set).
- To clear:
 - Press CLEAR to reset timer and memory to 00:00:00.
 - Or press HOUR + CLEAR, MINUTE + CLEAR, or SECOND + CLEAR to reset digits individually.

Start/Stop Countdown

- After time setting is ready, press START/STOP to begin countdown (1-sec resolution).
- Press again to stop/resume.
- When timer reaches 0:00:00:
 - Buzzer sounds
 - Corresponding indicator flashes slower.
 - Automatic count-up begins.
 - If multiple timers reach 0:00:00, all indicators flash slower, but only the displayed timer sounds.
- Press any button to stop alarm and count-up.

Memory Recall

- Press START/STOP to recall previous timer setting.
- Press again to start with recalled setting.

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Stopwatch Mode (Count-Up)

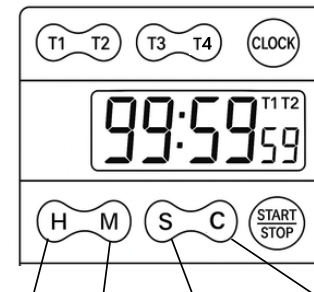
- In timer mode, press CLEAR to reset to 00:00:00.
- Press START/STOP to start/stop count-up (1-sec resolution).
- Max time: 99:59:59 → then auto-restarts from 00:00:00.

Placement Options

- Clip: Attach to board or pocket.
- Magnet: Stick to metal surfaces.
- Stand: Flip out rear metal stand for tabletop use.

Battery Replacement

- Turn timer over. Use a coin to open battery cover.
- Replace with 1 × 3V CR2032 (positive '+' facing up).
- Dispose of used battery according to local regulations.

HOUR MINUTE SECOND CLEAR

Distributed by Ultident Scientific | ULTIDENT.COM

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